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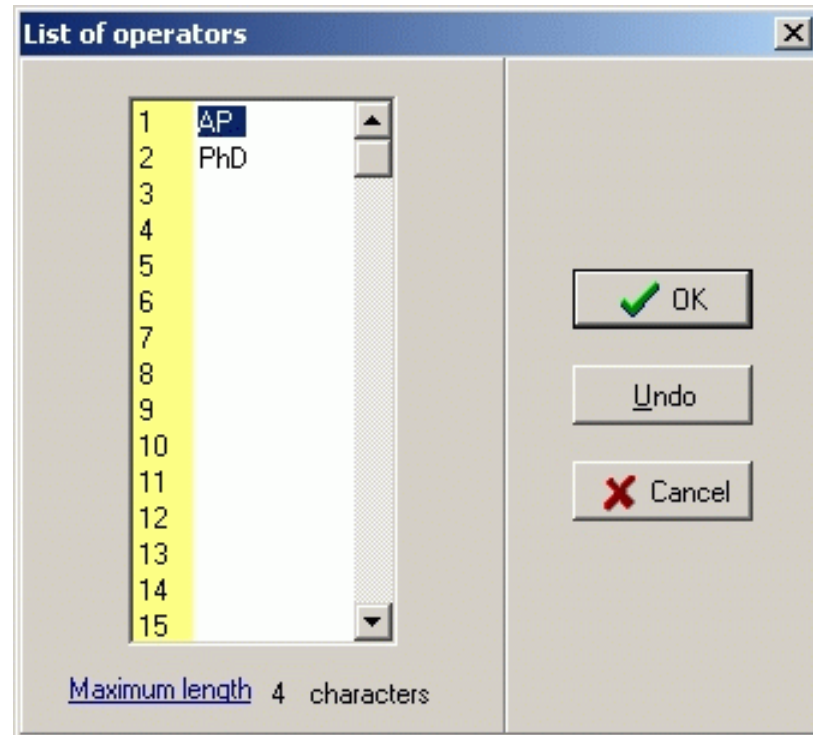
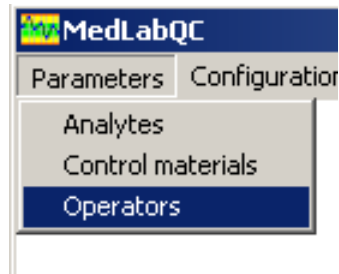
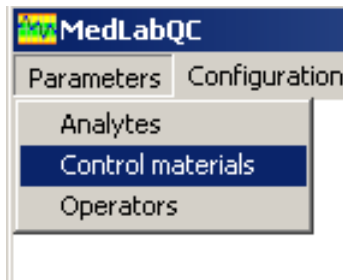
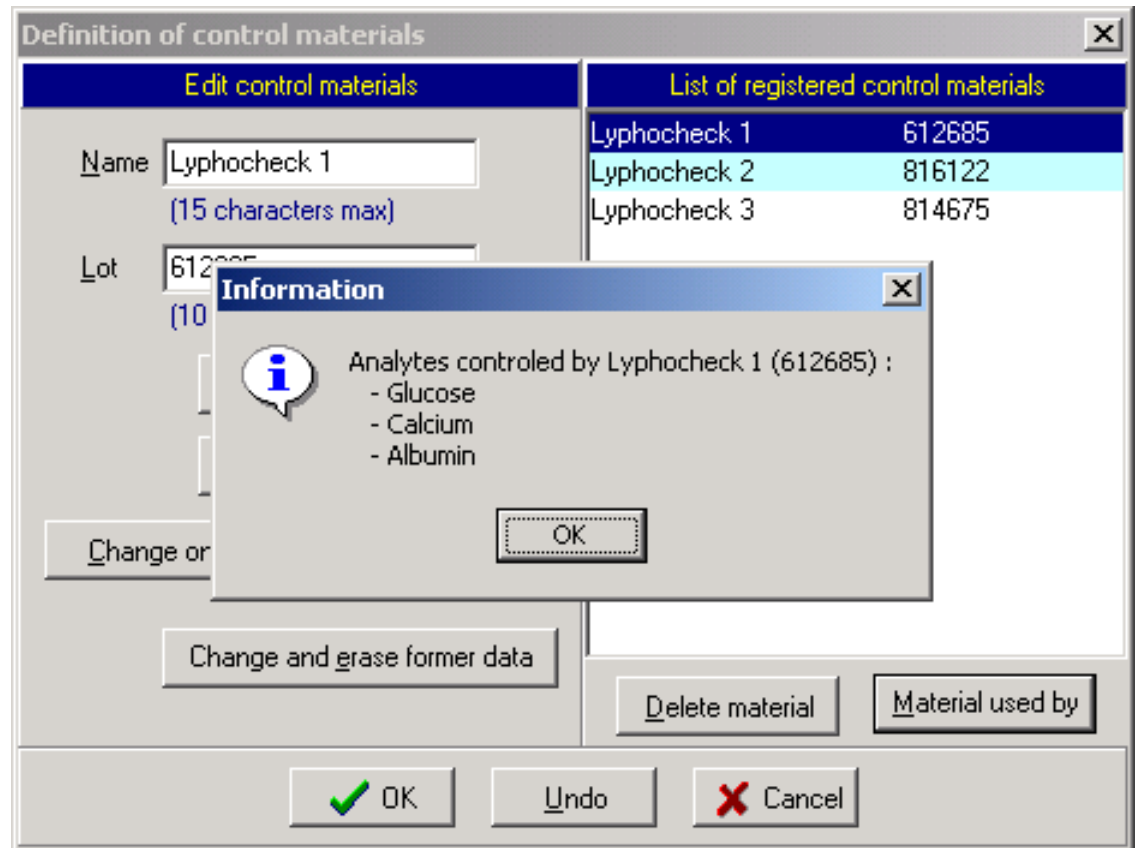
Creatinin (3 levels / $\mu\text{mol/l}$ / 0 decimals)

Iron (2 levels / $\mu\text{mol/l}$ / 1 decimal)

**→ Don't forget to click on "Add to list"
before clicking on "OK"**


ANALYTES BY CONTROL MATERIAL

Click on a control material and then click on « Material used by »

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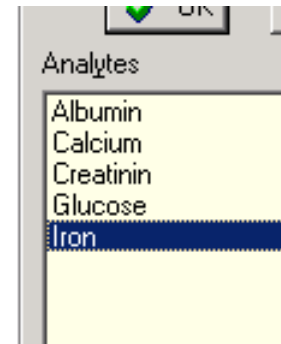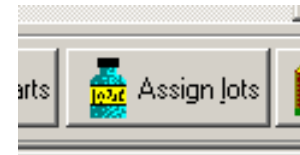
Validation

Change only the assignment of materials

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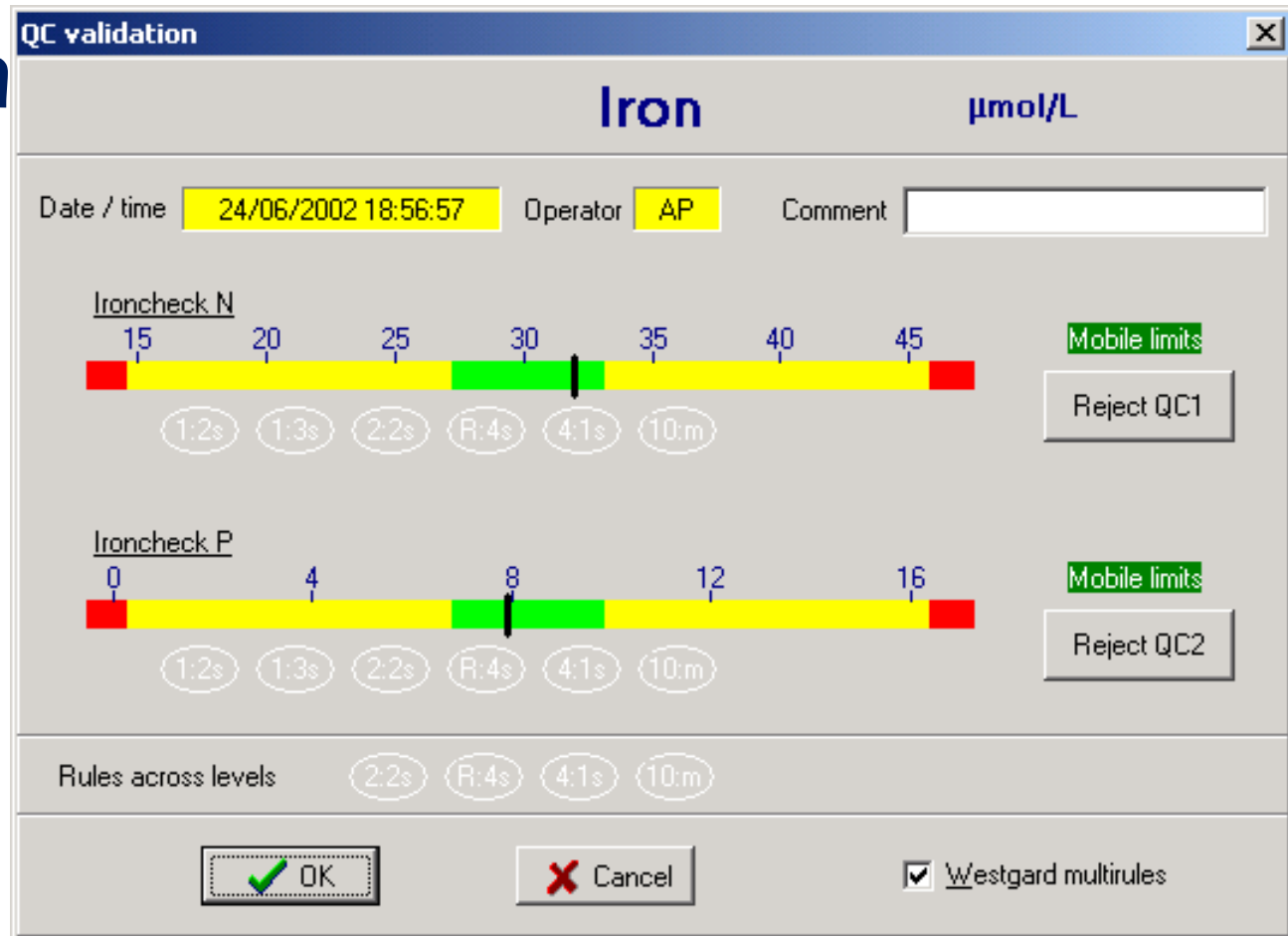
Westgard multirules

WESTGARD RULES (1/2)

1:2s = one observation in the warning zone

1:3s = one observation in the rejection zone

2:2s = 2 consecutive control observations in the warning zone, on the same side of the mean

R:4s = 2 consecutive control observations in the warning zone, on opposite sides of the mean

WESTGARD RULES (2/2)

4:1s = 4 consecutive control observations at a distance from the mean exceeding half the warning deviation

10m = 10 consecutive observations on the same side of the mean

→ There are other rules if you have more than one control level, named “rule across levels” for example:

R4s = 2 observations simultaneously in the warning zone, on opposite sides of the mean

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Locking control limits

From: 30/04/2002 to 15/05/2002

Partial statistics for the interval

#	m	CV	N
C1	2,377	4,6	16
C2	2,592	3,7	16

Lyphocheck 1
Locked from 30/04/2002 to 15/05/2002

Lyphocheck 2
Locked from 30/04/2002 to 15/05/2002

Fix →

Fixing control limits

E. coli

Target value: mm

Deviations allowed before
warning: action:

P. aeruginosa

Target value: mm

Deviations allowed before
warning: action:

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The screenshot shows the MedLabQC software interface. The 'Special functions' menu is open, with 'Lock limits' selected. The main window displays two control charts for Glucose. The top chart has a mean of 101.33 mg/dl and a CV of 1.5%. The bottom chart has a mean of 201.33 mg/dl and a CV of 1.6%. Both charts show data points for January 12, 2012, with values around 100 and 200 respectively. The right panel shows the 'Glucose' section with 'Locking control limits' and 'Mobile limits' options. The 'Mobile limits' section includes a table for partial statistics for the interval.

#	m	CV	N
C1			0
C2			0

Mobile limits
m = 101.33 mg/dl
CV = 1.5 %
N = 3

Mobile limits
m = 201.33 mg/dl
CV = 1.6 %
N = 3

locking control limits
From [] to []

Partial statistics for the interval

precinorm n
Mobile limits
Locked Mobile

precinorm p
Mobile limits
Locked Mobile

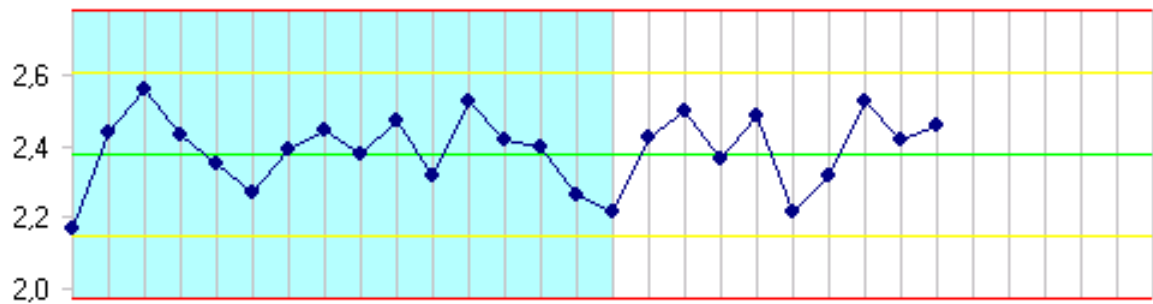
OK Undo Cancel

Data table Daily QC Control limits Print charts Assign jobs Archives Backup Queue : 2 Exit

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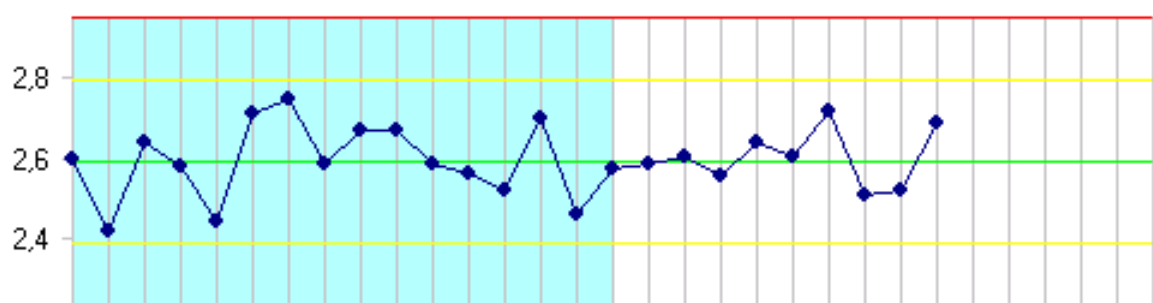
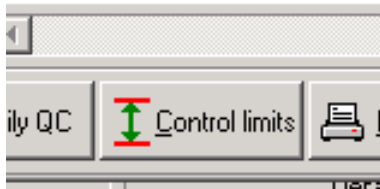
MedLabQC

Parameters Configuration File Clipboard **Special functions** ?

- Lock limits
- Fix limits**
- Erase/archive levels
- Random generator
- Clear queue

Mobile limits
m = 101.33 mg/dl
CV = 1.5 %
N = 3

Mobile limits
m = 201.33 mg/dl
CV = 1.6 %
N = 3

2 2 2
Jan
12

Glucose

Fixing control limits

precinorm n
Target value mg/dl
Deviations allowed before
warning action

precinorm p
Target value mg/dl
Deviations allowed before
warning action

OK

Data table Daily QC Control limits Print charts Assign lots Archives Backup Queue : 2 Exit